

Science.  
Applied to Life.™

3M Food Safety

# Simplicity is not the start, it's the result.

Get more than just results with  
3M™ Allergen Testing.



# The benefits are clear.

## 3M™ Allergen Protein Rapid Tests

### Lateral flow devices for specific allergen detection

Streamline your process and simplify your allergen testing with 3M™ Allergen Protein Rapid Tests – qualitative immunochromatographic assays for rapid in-plant monitoring of specific food allergens. Designed for accurate detection of processed and unprocessed allergen proteins, these fast, easy tests can be used for clean-in-place (CIP) final rinse water, environmental swab samples, raw ingredients and finished food products. We've taken the hassle out of allergen testing. You see the results.

### Lateral Flow Rapid Kits

Allergen protein	Catalogue	Assay time	Limit of detection
Almond	L25ALM	11 min ± 1 min	2.0 ppm
Bovine Total Milk	L25MLK	11 min ± 1 min	3.0 ppm
Cashew	L25CHW	11 min ± 1 min	2.0 ppm
Coconut	L25COC	11 min ± 1 min	2.0 ppm
Egg White	L25EGG	11 min ± 1 min	0.5 ppm
Fish	L25FSH	11 min ± 1 min	1.0 ppm
Gluten	L25GLU	11 min ± 1 min	5.0 ppm
Hazelnut	L25HZL	11 min ± 1 min	2.0 ppm
Peanut	L25PNT	11 min ± 1 min	1.0 ppm
Pecan	L25PEC	11 min ± 1 min	3.0 ppm
Pistachio	L25PST	11 min ± 1 min	2.0 ppm
Soy	L25SOY	11 min ± 1 min	2.0 ppm
Walnut	L25WAL	11 min ± 1 min	2.0 ppm

All kits include 25 lateral flow devices.






### Hook line assurance.

Our signature hook line feature can alert you to a possible false negative due to high amounts of target protein in a sample. This feature can increase confidence in your allergen testing results.

### Positive result .

Allergen detected in sample.

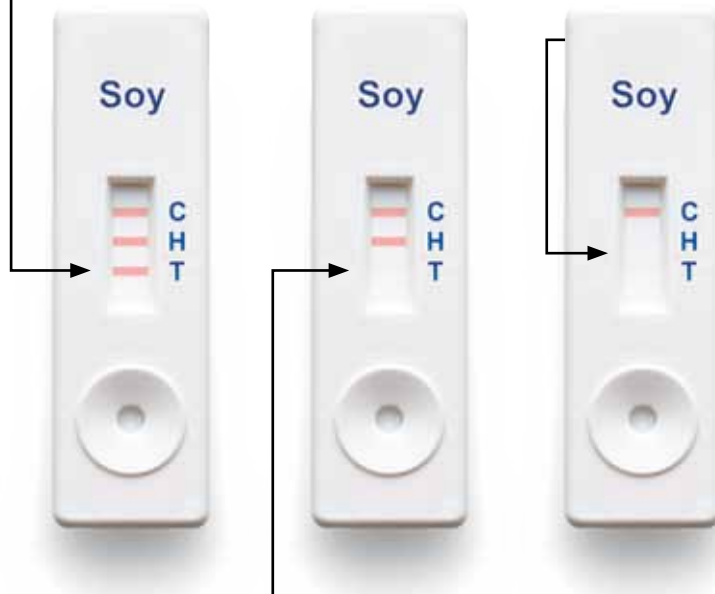

### Overloaded result.

If neither hook line nor test line appear, sample may be overloaded with too much allergen protein and should be diluted and re-tested.

### Negative result.

No allergen detected in sample.



info@hygiene-diagnostics.se • www.hygiene-diagnostics.se